

DATASHEET

Gelsolin (plasma isoform)

From porcine plasma
(>90% purity)

Cat. #: 8324-01
Quantity: 50µg

For Use in Research Only.
Not for Use in Diagnostic Processes.

Product Description

Plasma gelsolin has a molecular mass of ~80kD and consists of six homologous subdomains (S1-S6). Gelsolin is a Ca⁺⁺-dependent lipid binding actin regulatory protein, possessing three different actin modulating activities: severing of F-actin, nucleation of actin polymerization and capping of filaments barbed end. Gelsolin has three actin binding sites, two high affinity Ca⁺⁺ binding sites and two binding sites for PIP₂.

The protein concentration of gelsolin was determined by Biuret and the purity by scanning densitometry from Coomassie G-250 stained SDS-Gels. Lyophilized gelsolin (50µg) contains 10mM imidazole pH 7.0, 0.2mM DTT, 0.2mM EGTA, 2mM NaN₃ and 1% disaccharides, when reconstituted with 50µl H₂O.

The fragmentation activity of gelsolin was determined by Ostwald viscometry at 25°C. The specific viscosity of F-actin (1mg/ml) is reduced by ~65% after addition of gelsolin at a molar ratio of 1:100 (GS:A) after 30min. For this purpose the gelsolin stock solution was used as described below containing 1% disaccharides. F-actin containing 100mM KCl, 10mM Imidazole pH 7.4, 1mM ATP was mixed with 0.2mM CaCl₂ (final concentration) prior to the addition of gelsolin.

Product Handling

Preparation of a working stock

Add 50µl H₂O to the tube with gelsolin to obtain a 1mg/ml stock. Allow the protein to rehydrate for 2min and use a pipette to dissolve. Depending on the purpose this stock solution is ready-to-use. For critical assays we recommend buffer exchange against 10mM imidazole pH 7.0, 0.2mM DTT, 0.2mM EGTA and 2mM NaN₃.

Product Storage and Stability

Store the product at -70° upon arrival, where it will be stable for at least 6 months. Once dissolved, gelsolin is kept on ice and should be used within 5 days. Avoid refreezing

HYPERMOL EK WARRANTS THAT ITS PRODUCT ARE CONFORM TO THE INFORMATION CONTAINED IN THIS PUBLICATION. THE PURCHASERS MUST DETERMINE THE SUITABILITY OF THE PRODUCT(S) FOR THEIR PARTICULAR USE. ADDITIONAL TERMS AND CONDITIONS MAY APPLY (SEE REVERSE SIDE OF THE INVOICE). HYPERMOL BRANDS ARE SOLD EXCLUSIVELY BY HYPERMOL EK, GERMANY AND AUTHORIZED DISTRIBUTORS.

1 of 1

For product inquiries please contact:

cusserv@hypermol.com
Fon: +49 (0)521 9876228
Fax: +49 (0)521 9876231
www.hypermol.com