

## DATASHEET

### Biotin Actin

From rabbit skeletal muscle  
( $\alpha$ -actin)

For Use in Research Only.  
Not for Use in Diagnostic Processes.

Quantity: 2x100 $\mu$ g  
Cat.#: 8109-01

#### Product Description

G-actin from rabbit skeletal muscle is a single chain polypeptide with a molecular weight of 42kD consisting of 375 amino acids. Biotin-Actin is chemically modified G-actin by esterification with Biotin. Choosing a biotinylation using a  $\sim$ 20 Å spacer, renders the Biotin group to be readily available for antibody detection. Biotin-Actin is >99% pure according to scanning densitometry from Coomassie Blue G-250 stained PAA-Gels.

Biotin Actin is validated to polymerize in the presence of 100mM KCl, 2mM MgCl<sub>2</sub>. Immunoblotting using ExtrAvidin® alkaline phosphatase (Sigma, Cat. #: E2636) 1:10,000 with JetStain (Cat. #: 6022) detected 1ng of Biotin Actin.

Biotin-Actin is supplied as a lyophilized powder containing 2mM Tris-Cl pH 8.2, 0.4mM ATP and 0.1mM DTT, 1mM NaN<sub>3</sub>, 0.2mM CaCl<sub>2</sub> and 0.5% disaccharides when reconstituted with H<sub>2</sub>O to a 1mg/ml (23.8 $\mu$ M) solution.

For dilution of G-actin or exchange of ligand buffer into actin compatible buffer, MonoMix (Cat no.: 5100-0\*) may be used. In the absence of nucleators, the polymerization of G-actin can be initiated by PolyMix (Cat no.: 5000-0\*).

#### Product Handling

##### Preparation of a working stock

Example: Add 100 $\mu$ l of H<sub>2</sub>O to each tube containing 100 $\mu$ g Biotin-Actin to obtain a working stock of 1mg/ml (23.8 $\mu$ M). Allow the G-actin solution to rehydrate for 5 minutes at room temperature and vortex mildly or and dissolve by pipetting up and down. Leave at RT for another 5 min for rehydration. The final concentration of Biotin G-actin should not exceed 3mg/ml. For standard applications the Biotin-Actin working stock is now ready to use.

For critical assays the working stock should be applied to a gel filtration column or dialyzed (100vol. of MonoMix, O/N) followed by high-speed centrifugation. For less critical assays the actin working stock may be pre-spun (15.000rpm, 10min).

HYPERMOL EK WARRANTS THAT ITS PRODUCTS ARE CONFORM TO THE INFORMATION CONTAINED IN THIS PUBLICATION. THE PURCHASERS MUST DETERMINE THE SUITABILITY OF THE PRODUCT(S) FOR THEIR PARTICULAR USE. ADDITIONAL TERMS AND CONDITIONS MAY APPLY (SEE REVERSE SIDE OF THE INVOICE). HYPERMOL BRANDS ARE SOLD EXCLUSIVELY BY HYPERMOL EK, GERMANY AND AUTHORIZED DISTRIBUTORS.

1 of 1

*For product inquiries please contact:*

cusserv@hypermol.com  
Fon: +49 (0)521 9876228  
Fax: +49 (0)521 9876231  
www.hypermol.com